


How to adjust an Eppendorf Research Plus Fixed Pipette

(Figure 1):

1. Remove the ejector button (A) by pushing firmly upward.
2. Remove the cover panel with the color band (B) by pushing in on the left side and then push up from the bottom.
3. Remove the ring (C) and plunger button (D) by pulling firmly upward.

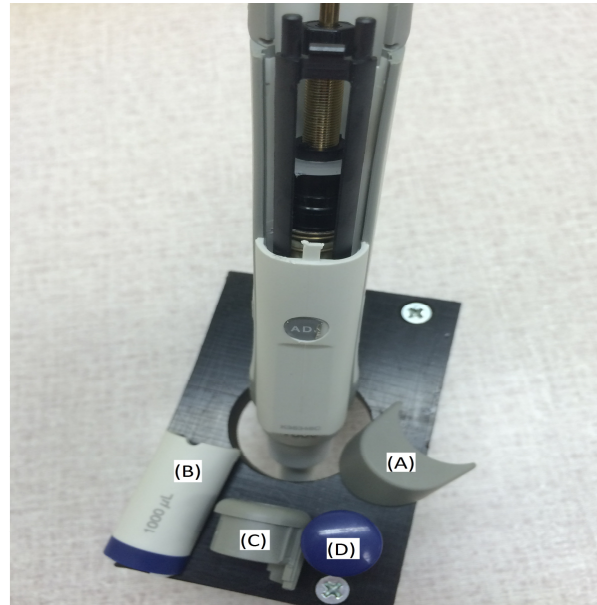

(Figure 1)

(Figure 2):


4. Locate the adjustment screw inside the unit.



(Figure 2)

(Figure 3):

5. Place adjustment tool over the adjustment screw and turn counter-clockwise to increase, and clockwise to decrease.
6. Remove adjustment tool, and replace all parts that were previously removed.

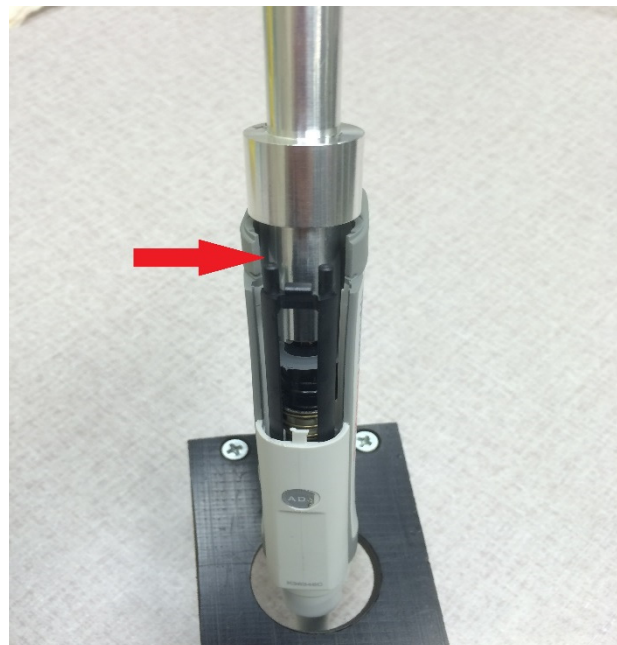

(Figure 3)