

macroman™

ADAPTOR FOR PIPETMAN® DIAMOND TIPS D5000 AND D10ML


MACROMAN™ can be used with PIPETMAN® DIAMOND tips D5000 and D10mL to offer a sterile and economical solution to transfer liquid.

This package includes:

- Tip adaptor
- Sample pack of D5000
- Sample pack of D10mL
- Bag of 10 filters

How to use it:

- Gently push the tip adaptor into the pipette holder and ensure it fits securely.
- Equip the adaptor with a polypropylene filter that prevents contamination.
- Fit either a D5000 or D10mL tip.

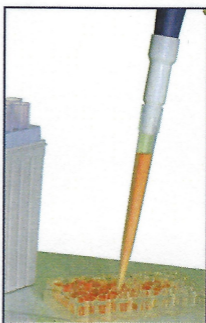

ORDERING INFORMATION

TYPE	PART NUMBER
EasyPack D5000 (5 x 200 tips)	F123969
Tipack D5000 (12 x 50 tips)	F161370
EasyPack D10mL (4 x 50 tips)	F161210
Tipack D10mL (8 x 30 tips)	F161230
Bag of 10 filters	F161280
Bag of 100 filters	F161241

 **GILSON®**