# seripettor® · seripettor® pro

## Mode d'emploi

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## Operating Manual

Please read carefully before use, and follow all operating and safety instructions!

BrandTech Scientific, Inc. · 11 Bokum Road · Essex, CT 06426-1506 · USA  
Phone: (860) 767-2562 · Fax: (860) 767-2563 · E-Mail: [mail@brandtech.com](mailto:mail@brandtech.com) · Internet: [www.brandtech.com](http://www.brandtech.com)

<b>Inhalt</b>		<b>Contents</b>		<b>Sommaire</b>		<b>Contenido</b>	
	<b>Seite</b>		<b>Page</b>		<b>Page</b>		<b>Página</b>
<b>Sicherheitsbestimmungen</b>	4	<b>Safety Instructions</b>	4	<b>Règles de sécurité</b>	4	<b>Normas de seguridad</b>	4
<b>Funktion</b>	6	<b>Application</b>	6	<b>Fonction</b>	6	<b>Función</b>	6
<b>Materialien</b>	7	<b>Materials</b>	7	<b>Matériaux</b>	7	<b>Materiales</b>	7
<b>Hinweis zur Geräteauswahl</b>	7	<b>Selecting the right dispenser</b>	7	<b>Sélectionner un distributeur</b>	7	<b>Información sobre la selección de instrumento</b>	7
<b>Einsatzausschlüsse</b>	8, 9	<b>Operating Exclusions</b>	8, 9	<b>Interdictions d'emploi</b>	8, 9	<b>Excepciones de uso</b>	8, 9
<b>Einzelteile, Montage, Demontage</b>	10, 11	<b>Components, Assembling, Dismantling</b>	10, 11	<b>Pièces détachées, montage, démontage</b>	10, 11	<b>Componentes, montaje, desmontaje</b>	10, 11
<b>Gerät zum Einsatz vorbereiten</b>	12	<b>Preparing instrument for use</b>	12	<b>Préparation de l'appareil pour l'emploi</b>	12	<b>Preparación del aparato para su uso</b>	12
Gerät entlüften und Volumen einstellen	13	Priming the instrument and setting the volume	13	Désaéragé et réglage du volume	13	Purgar el aire y ajustar el volumen	13
<b>Dosieren</b>	14	<b>Dispensing</b>	14	<b>Dosage</b>	14	<b>Dosificación</b>	14
<b>Wechsel der Dosiereinheit</b>	16	<b>Replacing the dispensing cartridge</b>	16	<b>Changement de l'unité de dosage</b>	16	<b>Cambio de la unidad de dosificación</b>	16
<b>Reinigung</b>	17	<b>Cleaning</b>	17	<b>Nettoyage</b>	17	<b>Limpieza</b>	17
Standardreinigung	18	Normal cleaning procedure	18	Nettoyage standard	18	Limpieza estandard	18
Ventile reinigen	19	Cleaning valves	19	Nettoyage des soupapes	19	Limpieza válvulas	19
<b>Störung – was tun?</b>	20	<b>Troubleshooting</b>	21	<b>Dérangement – que faire?</b>	22	<b>¿Qué hacer en caso de avería?</b>	23
<b>Technische Daten/ Bestelldaten</b>	24	<b>Technical Data/ Ordering Data</b>	24	<b>Données techniques/ Données de commande</b>	24	<b>Datos técnicos/ Referencias</b>	24
<b>Lieferumfang</b>	24	<b>Items supplied</b>	24	<b>Emballage standard</b>	24	<b>Alcance del suministro</b>	24
<b>Zubehör und Ersatzteile</b>	25, 26	<b>Accessories and Spare Parts</b>	25, 26	<b>Accessoires et pièces de rechange</b>	25, 26	<b>Accesorios y piezas de recambio</b>	25, 26
<b>Zur Reparatur einsenden</b>	27	<b>Return for Repair</b>	27	<b>Retour pour réparation</b>	27	<b>Envíos para reparación</b>	27
<b>Mängelhaftung</b>	28	<b>Warranty</b>	28	<b>Garantie</b>	28	<b>Garantía</b>	28
<b>Leitfaden zur Geräteauswahl</b>	29	<b>Dispenser selection chart</b>	30	<b>Table de sélection de distributeur</b>	31	<b>Tabla de selección de dispensadores</b>	32

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Read carefully before use!

## Safety Instructions

This instrument may be involved with hazardous materials, operations, and equipment.

This manual does not purport to address all of the safety risks associated with its use. It is the responsibility of whomever uses this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Follow the general instructions for hazard prevention and safety regulations, e.g., wear protective clothing, eye protection and gloves.
2. Observe the reagent manufacturer's information.
3. Every user must be acquainted with this operating manual before operation.
4. Use the instrument only for dispensing liquids, with strict regard to the defined Operating Exclusions (► pages 8, 9).  
If in doubt, contact the manufacturer or supplier.

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## Application

This instrument is designed for dispensing liquids, observing the following limits:

- vapor pressure up to 500 mbar
  - density up to 2.2 g/cm<sup>3</sup>
  - temperature 15 to 40 °C (seripettor® agar cultures up to 60 °C)
  - viscosity
    - 10 ml instrument: 150 mm<sup>2</sup>/s
    - 25 ml instrument: 75 mm<sup>2</sup>/s
- (dynamic viscosity [mPas]  
= kinematic viscosity [mm<sup>2</sup>/s]  
x density [g/cm<sup>3</sup>])

Valve block, filling tube, discharge tube and valves are autoclavable at 121 °C.

The dispensing cartridge (E) and pump assembly (A/D) are not autoclavable.

### Warning!

If there is a sign of a potential malfunction (e.g., piston difficult to move) never use force. Immediately stop dispensing and follow cleaning instructions (► page 17) or contact the manufacturer.

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## Materials

When the instrument is correctly used, the dispensed liquid comes into contact with only the following materials:

- seripettor®: PP, PE, EPDM
- seripettor® *pro*: PP, PE, borosilicate glass, Al<sub>2</sub>O<sub>3</sub>, ETFE, PFA, FEP, PTFE, Pt-Ir

Please see our catalog for a listing of chemical compatibility with components of this instrument. For further information please visit our internet site [www.brand.de](http://www.brand.de) or contact BRAND.

## Selecting the right dispenser

See "Dispenser selection chart" (► pages 29 - 32) for guidelines on selecting the right dispenser.

### Note:

Consult the following chapters for operating exclusions with your application (► pages 8, 9).

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### ■ seripettor®

#### Never use this instrument with

- liquids attacking PP, PE or EPDM
- non-polar solvents like hydrocarbons and halogenated hydrocarbons
- concentrated or oxidizing acids
- explosive liquids (e.g., carbon disulfide)

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### ■ seripettor® pro

#### Never use this instrument for

- liquids attacking PP, PE,  $Al_2O_3$ , ETFE, FEP, PFA and PTFE (e.g., dissolved sodium azide)
- liquids attacking borosilicate glass (e.g., hydrofluoric acid)
- liquids which are decomposed catalytically by platinum-iridium (e.g.,  $H_2O_2$ )
- non-polar solvents like hydrocarbons and halogenated hydrocarbons
- concentrated or oxidizing acids (excluding HCl)
- explosive liquids (e.g., carbon disulfide)
- suspensions (e.g., of charcoal) as solid particles may clog or damage the instrument

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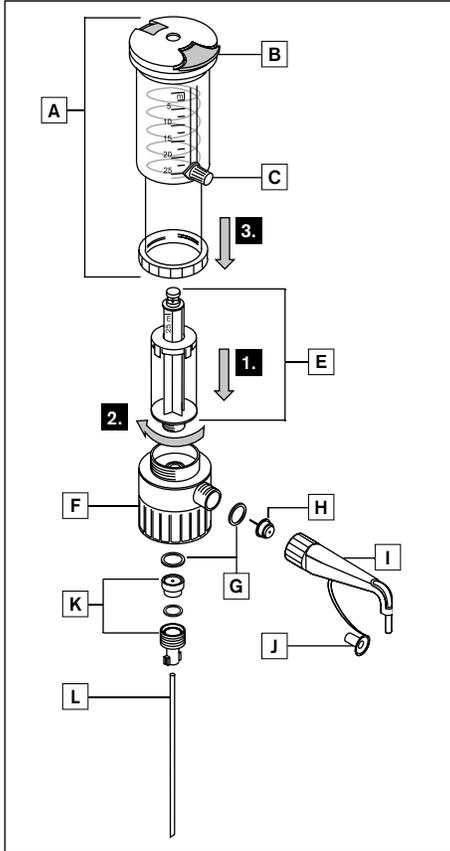
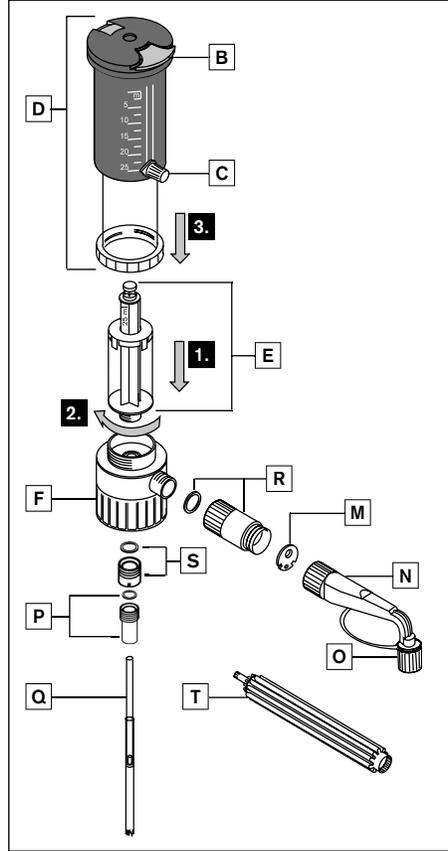
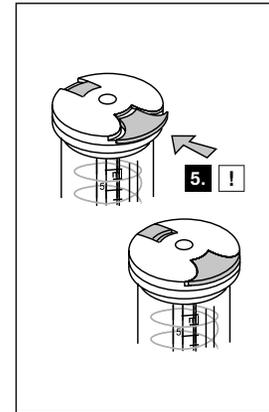
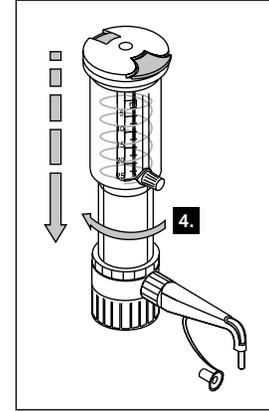
## ■ seripettor®

- A Pump assembly [PC]
- B Piston lock
- C Volume adjustment
- E\* Dispensing cartridge  
[piston: PP, cylinder: PE]
- F\* Valve block [PP]
- G\* Seals [EPDM]
- H\* Discharge valve [PP, EPDM]
- I\* Discharge tube [PP]
- J\* Stopper cap [PP]
- K\* Filling valve [PP] with O-ring  
[EPDM]
- L\* Filling tube [PP]

## ■ seripettor® pro

- D Pump assembly [PPO, PE]
- B Piston lock
- C Volume adjustment
- E\* Dispensing cartridge  
[piston: PP, cylinder: PE]
- F\* Valve block [PP]
- M Seal [PTFE]
- N\* Discharge tube with integrated valve  
[ETFE, FEP, PFA, Boro 3.3, Al<sub>2</sub>O<sub>3</sub>,  
Pt-Ir]
- O\* Stopper cap for discharge tube [PP]
- P\* Filling valve [ETFE, Boro 3.3, Al<sub>2</sub>O<sub>3</sub>]  
with seal [PTFE]
- Q\* Telescopic filling tube [FEP, PTFE]
- R\* Adapter for discharge tube [PP]  
with seal [PTFE]
- S\* Adapter for filling valve [PP]  
with seal [PTFE]
- T Mounting tool

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## Preparing instrument for use

### **Warning!**

Wear protective clothing, eye protection and gloves. Follow all Safety Instructions and observe the Operating Exclusions (► pages 4 - 5, 8 - 9).

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### **Note:**

Follow all instructions on pages 10, 11 for proper assembly.

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## Priming the instrument and setting the volume

### Attention!

To avoid splashes when priming, place the tube opening against the inner wall of a suitable vessel. For small bottles use a bottle stand to prevent tipping over.

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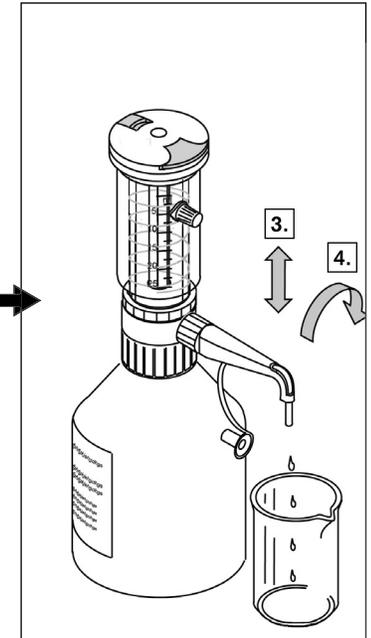
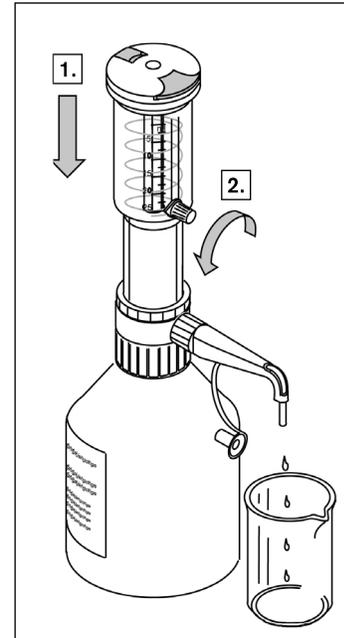
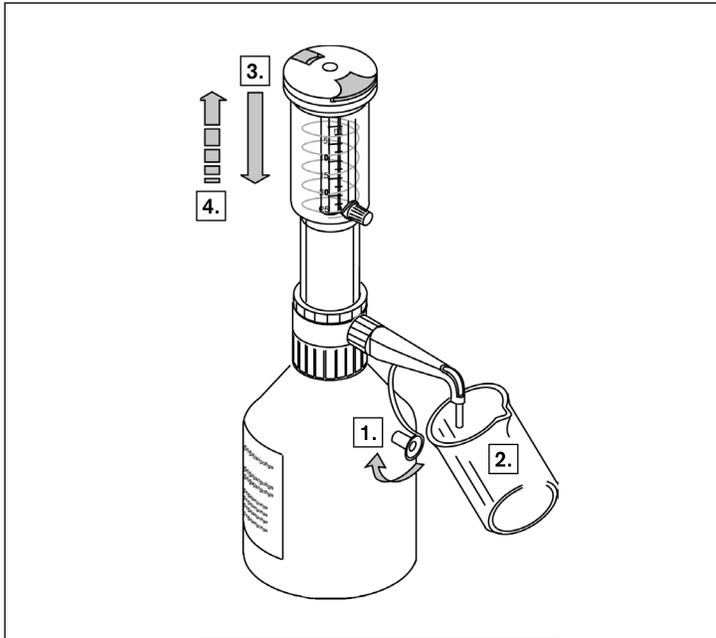
## Dispensing

### Attention!

Follow all Safety Instructions and  
observe the Operating Exclusions  
(▶ pages 4 - 9).

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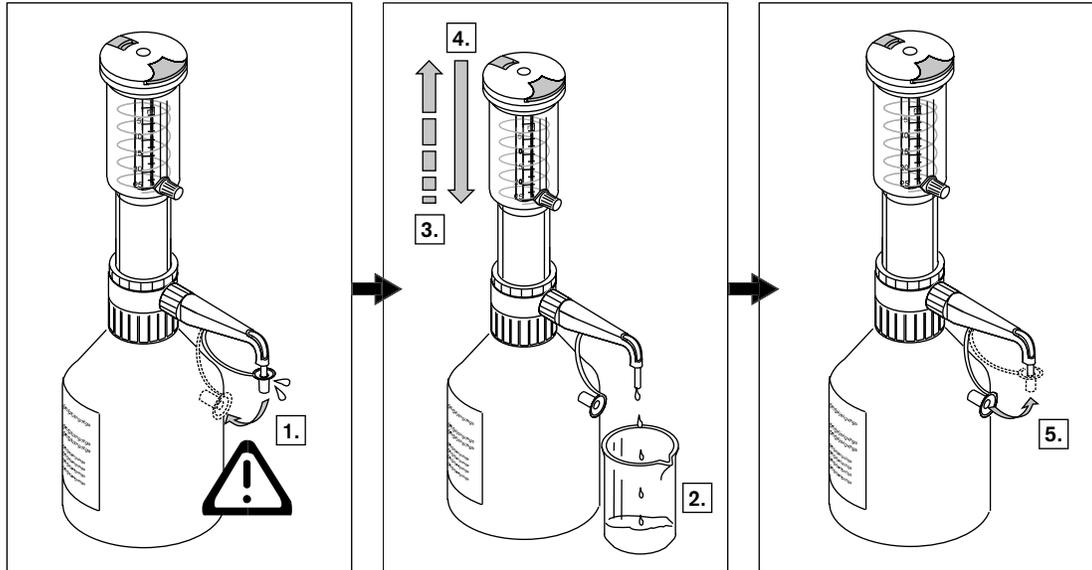
## Serial dispensing with flexible discharge tube (optional)

### Attention!

Each time you are going to use the tubing, examine it carefully for any damage.

To dispense aggressive liquids, e.g. concentrated acids and alkalis, you should take safety measures in addition to the normal precautions of wearing protective clothing, eye protection and protective gloves. We recommend that you use a protective shield.

When using the flexible discharge tube, the bottle must be supported using a bottle stand to help prevent the bottle from accidentally tipping over. To help avoid reagent splashing from the tube, always grip the tube firmly by the handle and replace in the holder after use.

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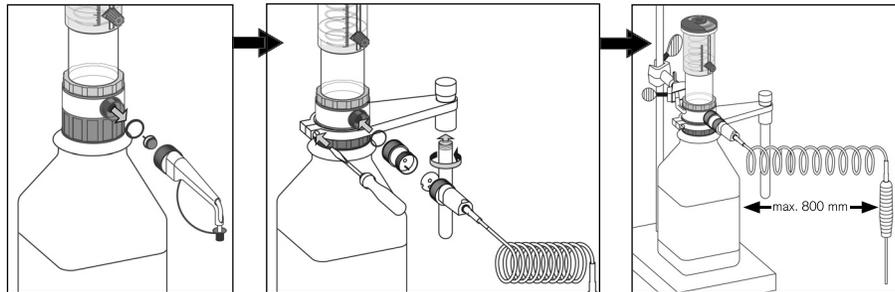
## Replacing the dispensing cartridge

### Attention!

The dispensing cartridge is a wearing part. Replace the cartridge when the piston is difficult to move or is not airtight.  
To prevent chemical injury, the dispensing cartridge has to be cleaned (► page 18) prior to replacement.

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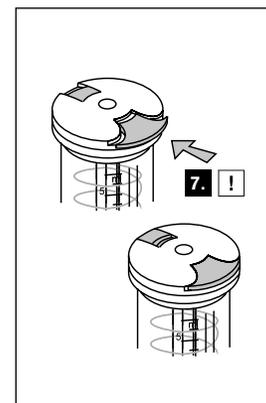
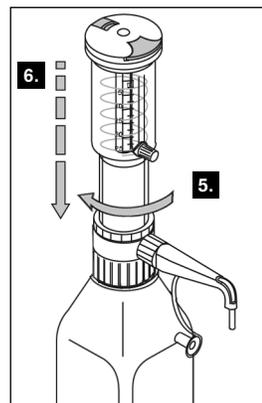
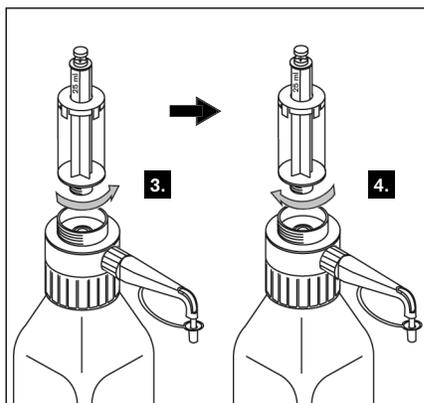
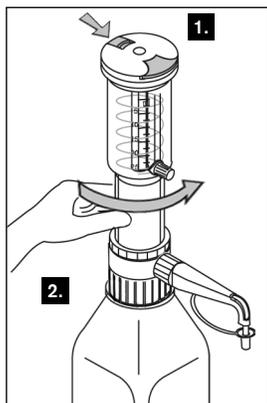
## Cleaning

The instrument must be cleaned in the following situations to assure correct operation:

- immediately, if the piston becomes difficult to move
- before changing the reagent
- prior to long term storage
- before replacing the valves
- before replacing the dispensing unit

Depending upon the frequency of use, the dispensing cartridge should be replaced regularly.

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### Warning!

Follow all Safety Instructions (► page 4, 5).

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#### Warning!

Carry the mounted instrument only as shown in the figure.

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#### Note:

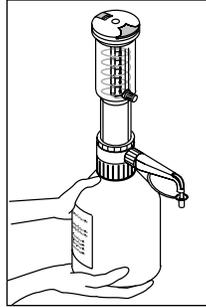
Clean only inside a suitable collecting basin.

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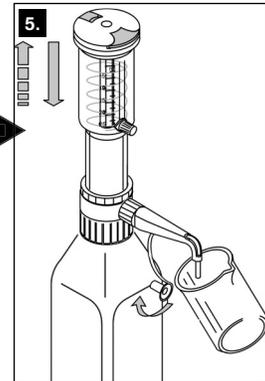
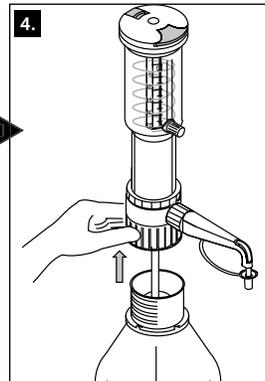
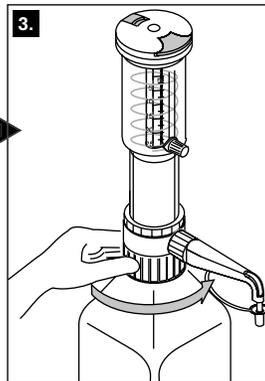
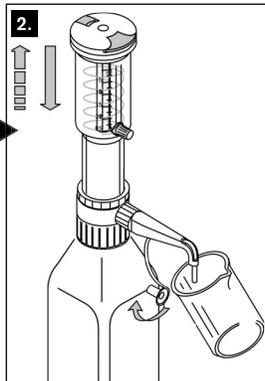
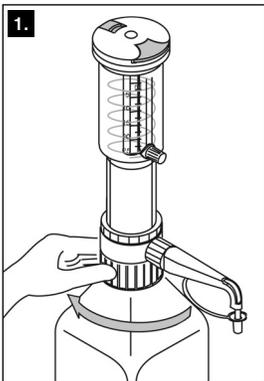
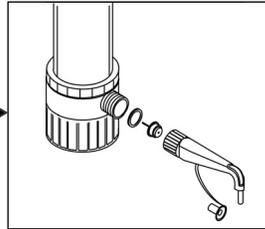
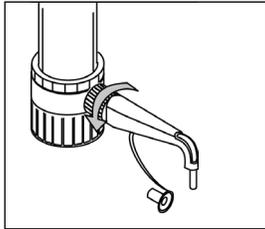
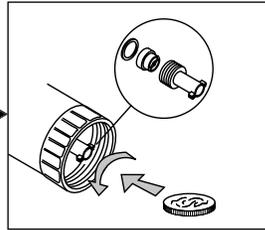
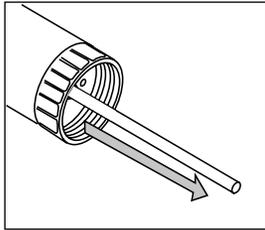
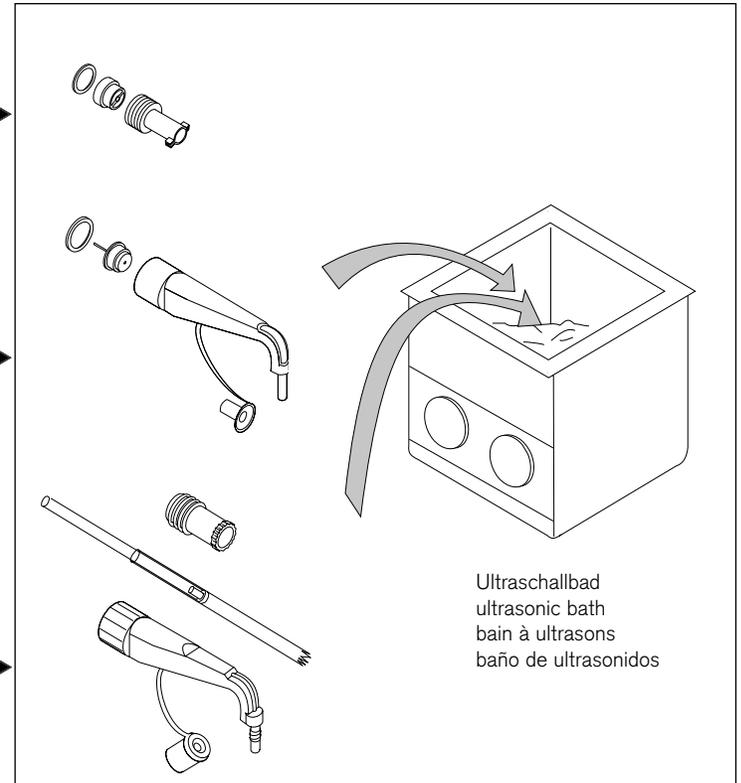
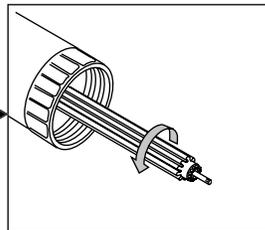
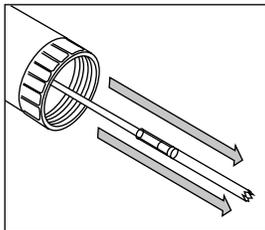
## Troubleshooting

Problem	Possible cause	Corrective action
<b>Piston is difficult to move</b>	<ul style="list-style-type: none"> <li>- Formation of crystals</li> <li>- Dispensing of incompatible liquid</li> </ul>	<ul style="list-style-type: none"> <li>- Replace dispensing cartridge.</li> <li>- Check for operating exclusions (▶ page 30).</li> </ul>
<b>Instrument does not draw in liquid</b>	<ul style="list-style-type: none"> <li>- Valve not correctly mounted</li> <li>- Wrong dispensing cartridge</li> <li>- Volume adjusted to minimum setting</li> <li>- Filling tube blocked or not mounted correctly</li> <li>- Sticking filling valve</li> <li>- Dispensing cartridge worn</li> </ul>	<ul style="list-style-type: none"> <li>- Mount valve correctly.</li> <li>- Mount compatible dispensing cartridge.</li> <li>- Set to required volume (▶ page 13).</li> <li>- Push the filling tube on firmly, cut off the end diagonally.</li> <li>- Clean filling valve (▶ page 19). If necessary, replace the valve and sealing washer (observe Operating Exclusions ▶ pages 8 - 9).</li> <li>- Replace dispensing cartridge (▶ page 16).</li> </ul>
<b>Instrument does not dispense liquid</b>	<ul style="list-style-type: none"> <li>- Sticking discharge valve</li> </ul>	<ul style="list-style-type: none"> <li>- Clean discharge valve. If necessary replace discharge valve.</li> </ul>
<b>Air bubbles in the instrument</b>	<ul style="list-style-type: none"> <li>- Not primed properly</li> <li>- Dispensing cartridge worn</li> <li>- Reagent with high vapor pressure has been drawn in too quickly</li> <li>- Air not expelled from the instrument</li> <li>- Filling tube not firmly connected or damaged</li> <li>- Valves not firmly connected or damaged</li> </ul>	<ul style="list-style-type: none"> <li>- Prime the instrument (▶ page 13).</li> <li>- Replace dispensing cartridge (▶ page 16).</li> <li>- Release piston slowly.</li> <li>- Purge the air (▶ page 13).</li> <li>- Follow normal cleaning procedure (▶ page 18). Push the filling tube on firmly. If necessary, cut off approx. 1 cm of tube at the upper end and re-connect it or replace filling tube.</li> <li>- Follow normal cleaning procedure (▶ page 18). Tighten the valves. If necessary, replace the valves and sealing washers.</li> </ul>
<b>Dispensed volume is too low</b>	<ul style="list-style-type: none"> <li>- Discharge tube not firmly connected or damaged</li> <li>- Filling tube not firmly connected or damaged</li> <li>- Valves not firmly connected or damaged</li> </ul>	<ul style="list-style-type: none"> <li>- Push the original manufacturer's discharge tube on firmly. Replace a deformed or damaged discharge tube.</li> <li>- Normal cleaning procedure (▶ page 18). Push the filling tube on firmly. If this does not suffice, cut off approx. 1 cm of the tube at the upper end and re-connect it or replace filling tube.</li> <li>- Clean the instrument (▶ page 18). Tighten the valves. If necessary, replace valves and sealing washers.</li> </ul>
<b>Liquid above piston seal</b>	<ul style="list-style-type: none"> <li>- Piston damaged</li> </ul>	<ul style="list-style-type: none"> <li>- Replace dispensing cartridge (▶ page 16).</li> </ul>

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\* The values of accuracy and coefficient of variation are final test values referring to the delivered volume, instrument and distilled water at equilibrium with ambient temperature (20 °C/60 °F) and smooth and steady operation.

A = Accuracy  
CV = Coefficient of variation

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## Warranty

As provided by law, any and all warranties are null and void if the product has been misused, modified or repaired by unauthorized personnel, if the defects are caused by negligence (non-observance of operating manual, lack of maintenance) or by normal wear and tear. Use only original manufacturer's accessory/spare parts.

For conditions and extent of warranty refer to our warranty card enclosed with the product.

Subject to technical modification without notice. Errors excepted.

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Medium	seripetto®	seripetto® <i>pro</i>
Formamide	+	+
Formic acid, 100%		+
Glycerol		+
Glycol (Ethylene glycol)	+	+
Glycolic acid, 50%	+	+
Hexanoic acid	+	+
Hexanol		+
Hydrolic acid	+	+
Hydrobromic acid		+
Hydrochloric acid, 37 %		+
Hydrogen peroxide, 35%	+	
Isamyl alcohol		+
Isobutanol	+	+
Isopropanol (2-Propanol)	+	+
Lactic acid	+	+
Methanol	+	+
Methyl benzoate		+
Methyl propyl ketone		+
Mineral oil (Engine oil)		+
Monochloroacetic acid		+
Nitric acid, 10%		+
Oxalic acid	+	+
Perchloric acid		+
Phenol		+
Phosphoric acid, 85%		+
Piperidine		+
Potassium chloride	+	+
Potassium dichromate	+	+
Potassium hydroxide	+	+
Potassium permanganate	+	+
Propionic acid	+	+
Propylene glycol (Propandiol)	+	+
Pyridine		+
Pyruvic acid	+	+
Salicylaldehyde		+
Salicylic acid	+	+
Silver acetate	+	+
Silver nitrate	+	+
Sodium acetate	+	+
Sodium chloride	+	+
Sodium dichromate	+	+
Sodium fluoride	+	+
Sodium hydroxide, 30%	+	+
Sodium hypochlorite	+	+
Sulfuric acid, 10%	+	+
Tartaric acid		+
Urea	+	+
Zinc chloride, 10%	+	+
Zinc sulfate, 10%	+	+

## seripetto® and seripetto® *pro* are not suitable for hydrofluoric acid (HF)

The above recommendations reflect testing completed prior to publication. Always follow instructions in the operating manual of the instrument as well as the reagent manufacturer's specifications. In addition to these chemicals, a variety of organic and inorganic saline solutions (e.g., biological buffers), biological detergents and media for cell culture can be dispersed. Should you require information on chemicals not listed, please feel free to contact BRAND. Status as of: 12/06

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